# UV LIGHT SOURCE

## USER MANUAL

Manual version: 1.0.A  
Product code: G2008  
Product Version: 1.0

# Contents

- 1. Overview ..... 3
- 2. EU Declaration of Conformity ..... 4
- 3. Safety..... 3
  - 3.1 Warning ..... 3
  - 3.2 Use of Equipment..... 3
  - 3.3 Servicing ..... 3
- 4. Unpacking..... 4
  - 4.1 Packing List..... 4
  - 4.2 Damage Inspection..... 5
- 5. Operation..... 5

# 1. Overview

Ossila's USB powered UV light source is a simple and flexible light source for steady-state spectroscopy applications. It can be used in conjunction with a spectrometer to perform fluorescence measurements.

## Key features

- UV emission (360 nm – 370 nm approx.)
- LED source for stable output
- USB powered
- Compact design

## 2. EU Declaration of Conformity

### We

**Company Name:** Ossila BV

**Postal Address:** Biopartner 3 building, Galileiweg 8

**Postcode:** 2333 BD Leiden

**Country:** The Netherlands

**Telephone number:** +31 (0)71 3322992

**Email Address:** info@ossila.com

**declare that the DoC is issued under our sole responsibility and belongs to the following product:**

**Product:** UV White Light Source (G2008A1)

**Serial number:** G2008A1- xxxx

### Object of declaration:

UV Light Source (G2008A1)

**The object of declaration described above is in conformity with the relevant Union harmonisation legislation:**

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## 3. Safety

### 3.1 Warning

The Ossila UV Light Source is classified as being in Risk Group 2 (Moderate Risk) under the IEC/EN 62471 (Photobiological Safety of Lamps and Lamp Systems) Standard for actinic UV skin & eye hazard. The maximum permissible exposure time is 2400 seconds. **Do not look directly into the source. Avoid skin and eye exposure.**

### 3.2 Use of Equipment

The Ossila UV Light is designed to be used as instructed. It is intended for use under the following conditions:

- Indoors in a laboratory environment (Pollution Degree 2).
- Altitudes up to 2000m.
- Temperatures of 5°C to 40°C; maximum relative humidity of 80% up to 31°C.

### 3.3 Servicing

If servicing is required, please return the unit to Ossila Ltd. The warranty will be invalidated if:

- Modification or service has taken place by anyone other than an Ossila engineer.
- The Unit has been subjected to chemical damage through improper use.
- The Unit has been operated outside the usage parameters stated in the user documentation associated with the Unit.
- The Unit has been rendered inoperable through accident, misuse, contamination, improper maintenance, modification, or other external causes.

## 4. Unpacking

### 4.1 Packing List

The standard items included with the Ossila UV Light are:

- Ossila UV Light.
- USB type-C cable.

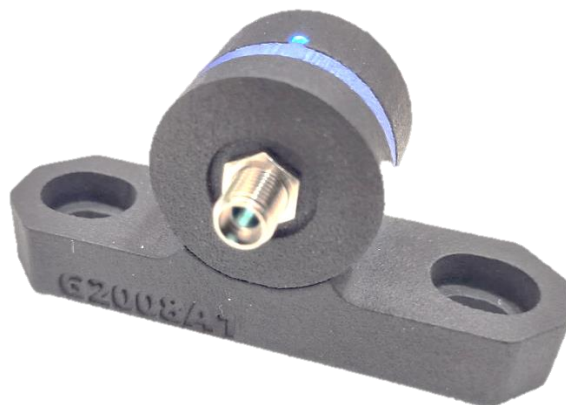
Figure 4.1. Ossila UV Light.

Figure 4.2. USB type-A to type-C cable.


## 4.2 Damage Inspection

Examine the components for evidence of shipping damage. If damage has occurred, please contact Ossila directly for further action.

## 5. Operation

The Ossila UV Light is powered by a USB type-C cable supplying 5 V at 100 mA. Any PC USB port or plug socket USB adapter can power the device. The source will turn on when the USB cable is plugged into the USB port on the rear of the device. A SMA905 optical fibre can be connected to the output port of the device. A blue indicator LED on the top of the unit will turn on while the system is powered. Do not look into the front of the device or expose any skin to the front of the device while the system is powered.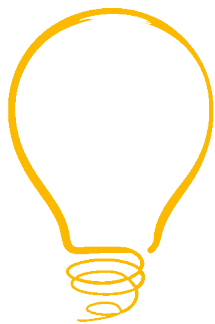

**INNOVATION  
AWARD 2017**

**IOP** Institute of Physics

THE QUEEN'S AWARDS  
FOR ENTERPRISE:  
INTERNATIONAL TRADE  
2018